AdvanBlock[™]-EIA Blocking Solution

Immunoassay blocking and antibody incubation solution

AdvanBlock[™]-EIA is a novel blocking solution, optimized to enhance specific antibody-antigen interactions for ELISA. This all-in-one blocking solution and antibody incubation buffer is designed to decrease non-specific binding caused by low quality antibodies and serum matrix effects. Provided as a convenient ready-to-use solution intended to directly replace other commonly used blocking buffers.

Advantages

- DECREASES BACKGROUND
- DECREASES WELL-TO-WELL VARIABILITY
- DECREASES NON-SPECIFIC BINDING
- READY-TO-USE SOLUTION



Figure 1. Decreased serum matrix effect with AdvanBlock[™]-EIA compared to 1% BSA/PBST. The ELISA plate was coated with Extractable Nuclear Antigens extract diluted in 1X EIA Coating Buffer. The plate was blocked with AdvanBlock[™]-EIA or 1% BSA-PBST before serial dilutions of Systemic Lupus Erythematosus plasma prepared in AdvanBlock[™]-EIA or 1% BSA-PBST were applied to the plate. The wells were washed with 1X AdvanWash[™] Buffer, then goat anti-human IgG/A/M-HRP diluted in AdvanBlock[™]-EIA or 1% BSA-PBST were added to the plate. The wells were washed and TMB substrate was added before stopping with sulfuric acid. Non-specific sample matrix effects and well-to-well variability that are observed with some patient samples are decreased with AdvanBlock[™]-EIA.

Ordering Information

| R-03728-E10 AdvanBlock [™] -EIA Blocking Solution 1 L | Catalog Number | Product | Size |
|--|----------------|--|------|
| | R-03728-E10 | AdvanBlock [™] -EIA Blocking Solution | 1 L |

Related Product

| K-16025-D10 | ELISABright™, sufficient for ten (10) 96-well ELISA plates | 100 ml |
|-------------|--|--------|
| R-03024-D50 | AdvanWash™ 10X Washing Solution | 500 ml |

Advansta Corporation

2140 Bering Drive | San Jose, CA 95131 Tel: 650.325.1980 | Fax: 650.325.1904 | Email: sales@advansta.com Product information: www.advansta.com/products/AdvanBlock-EIA Copyright © 2016 Advansta. All rights reserved. The Advansta logo is a registered trademark of the Company. AdvanBlock™, AdvanWash™ and ELISABright™ are trademarks of the Company. All other trademarks, service marks and tradenames appearing in this brochure are the property of their respective owners.



